# CAMCOLMS1

**1.3 MEGAPIXEL DIGITAL MICROSCOPE - 20-200X ZOOM**  
**1.3 MEGAPIXEL DIGITALE MICROSCOOP - 20-200X VERGROTING**  
**MICROSCOPE NUMÉRIQUE 1.3 MÉGAPIXELS - AGRANDISSEMENT 20-200X**  
**MICROSCOPIO DIGITAL 1.3 MEGAPÍXELES - AMPLIACIÓN 20-200X**  
**1.3 MEGAPIXEL DIGITALES MIKROSKOP - VERGRÖßERUNG 20-200X**  
**MICROSCÓPIO 1,3 MEGAPIXELS - 20-200X AMPLIAÇÃO**  
**MIKROSKOP CYFROWY - 1.3 MEGAPIKSELE - POWIĘKSZENIE 20 - 200X**

USER MANUAL	3
GEbruikersHANDLEIDING	9
MODE D'EMPLOI	15
MANUAL DEL USUARIO	21
BEDIENUNGSANLEITUNG	27
MANUAL DO UTILIZADOR	33
INSTRUKCJA OBSŁUGI	39

# USER MANUAL

## 1. Introduction

To all residents of the European Union

### Important environmental information about this product

This symbol on the device or the package indicates that disposal of the device after its lifecycle could harm the environment. Do not dispose of the unit (or batteries) as unsorted municipal waste; it should be taken to a specialized company for recycling. This device should be returned to your distributor or to a local recycling service. Respect the local environmental rules.

### If in doubt, contact your local waste disposal authorities.

Thank you for choosing Velleman! Please read the manual thoroughly before bringing this device into service. If the device was damaged in transit, don't install or use it and contact your dealer.

## 2. Safety Instructions

- The white LEDs which illuminate the microscope target field are very bright. Do not stare directly into these LEDs as it may damage your eyes.
- The clear plastic distance shell sometimes picks up dirt or toxic material from a microscopically observed surface. Be careful that this doesn't get in contact with the human skin. Always wash carefully or disinfect as needed.

## 3. General Guidelines

Refer to the **Velleman® Service and Quality Warranty** on the last pages of this manual.

- The microscope is not waterproof, so keep it dry.
- Do not use it in a humid place like bathrooms. A dry environment will maintain its life to maximum.
- Use the microscope only between -5°C and 50°C.
- Sudden temperature change may form dew inside the microscope. Put it inside a handbag or plastic bag to slowdown temperature changes.
- Do not point Microscope lens to the sun or strong light for a long time. Powerful light may hurt the light-sensitive electronics.
- Avoid touching the lens.
- Do not unplug the microscope from USB port when LEDs are on. This may cause information loss or circuit damage. Please always close the MicroCapture software and then unplug the microscope.

## 4. Features

- Ideal for checking PCBs, plants, insects, rocks, minerals...
- Adjustable focus and magnification
- IR LEDs with adjustable intensity

## 5. Overview

Refer to the illustrations on page 2 of this manual.

1	Snapshot button
2	Focus adjustment wheel

## 6. Operation

### 6.1 Installing the software

#### Microsoft Windows

To use the device, you need to install the following software from the included CD:

- The device driver
  - The Microcapture software, which allows using the microscope.
1. Insert the CD in your computer.  
The installation wizard will start automatically.  
If not, browse the CD through Windows Explorer and double-click the file **MicroCapture.exe**.
  2. The Microscope Installation window appears.

3. To install the device drive, click **USB Microscope Driver** and go through the steps of the wizard.
4. Next, click **MicroCapture software** to install the microscope software and go through the steps of the wizard.

- Restart if your computer asks you to.

### Apple Macintosh

- Insert the CD in your computer.
- Open the folder mac on the CD and copy MicroCapture onto your computer.

### 6.2 Starting the microscope


- Install the microscope on the stand.
- Use the included cable to connect the device to a USB port on your computer.
- Double click the Microcapture icon on your desktop to start the software.  
The software will start and display the image from the microscope.
- If necessary, use the wheel [2] on the device to adjust focus.

### 6.3 Capturing an image

- To capture an image, click the camera icon in the toolbar. Captured photos are saved automatically and image thumbnails will be available on the Photo tab on the right of the window for further editing.

### 6.4 Capturing video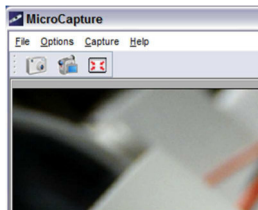

- To capture video, click the video camera icon in the toolbar.  
The icon will start flashing.
- To stop recording, click the icon again.  
Captured video is saved automatically and displayed on the Video tab on the right side of the window.

### Note

The most appropriate pixel setting for video capturing is 640x480 (see section Preview size below for settings).

## 6.5 Tasks for images

You can perform various tasks on the images you have captured.

- Click the Photo tab on the right side of the window.
- Right-click an image. Available tasks are:

Option	Function
Open	Opens the image in your default image programme
Save as	Saves the image with another name or format
Copy	Copies the image
Delete	Deletes the image

## 6.6 Tasks for video

You can perform various tasks on the videos you have captured.

- Click the Video tab on the right side of the window.  
The options in the context menu depend on your operating system.

## 6.7 Changing the default folder for captured images or video

You can change the default folder for your captured images or video.

- On the File menu, click Photos\_Directory (or Videos\_Directory for video).
- Select the folder you want and click OK.
- All new images will now be saved to that folder.

## 6.8 Software options

The Options menu contains the following options

Option	Function
<b>Preview size</b>	Set the preview size for pictures and video. The most appropriate pixel setting for video capturing is 640x480.
<b>Rotate</b>	Rotate the image clockwise or counter clockwise.

<b>Date/Time</b>	Show or hide the date and time on the preview
<b>Language</b>	Select another language for the software.
<b>Skin</b>	Select another skin for the software.
<b>FullScreen</b>	Set the microscope view to full screen. You can also use the Full Screen icon on the toolbar.
<b>Preferences</b>	Set preferences for pictures and video

## 7. Cleaning and maintenance

The clear plastic distance shell sometimes picks up dirt or toxic material from a microscopically observed surface. Be careful that this does not get in contact with the human skin. Always wash carefully or disinfect as needed.

## 8. Technical Specifications

image sensor	CMOS sensor
number of pixels	1.3 megapixel
effective pixels	1280 (H) x 1024 (V)
input interface	USB 2.0
magnification rate	20 x / 200 x
cable length	1.4 m
white balance	automatic
IR LEDs	8 LEDs (adjustable by control wheel)
iris aperture	AES
picture format	JPEG/BMP
picture	picture format: JPEG/BMP
	resolution: up to 1600 x 1200 pixels
video	video format: AVI
	resolution: up to 1600 x 1200 pixels
power supply	5 VDC via USB
system requirements	Windows® 2000/XP/Vista/Win7/MacOS®
dimensions	Ø 33 x 110 mm
weight	100 g

**Use this device with original accessories only. Velleman nv cannot be held responsible in the event of damage or injury resulting from (incorrect) use of this device.**

**For more info concerning this product and the latest version of this manual, please visit our website [www.velleman.eu](http://www.velleman.eu).**

**The information in this manual is subject to change without prior notice.**

All registered trademarks and trade names are properties of their respective owners and are used only for the clarification of the compatibility of our

products with the products of the different manufacturers. Windows, Windows XP, Windows 2000, Windows Vista, Windows Aero, Windows 7, Windows 8, Windows Mobile, Windows Server are registered trademarks of Microsoft Corporation in the United States and other countries. iPad, iPod, iPod touch, iPhone, Mac, iMac, MacBook, PowerBook, Power Mac, Mac OS are trademarks of Apple Inc., registered in the U.S. and other countries. Android is a trademark of Google Inc. The Trademark BlackBerry® is owned by Research In Motion Limited and is registered in the United States and may be pending or registered in other countries. Velleman NV is not endorsed, sponsored, affiliated with or otherwise authorized by Research In Motion Limited. Symbian is a registered trademark of Symbian Software Ltd, Symbian Ltd.

**© COPYRIGHT NOTICE**

**The copyright to this manual is owned by Velleman nv. All worldwide rights reserved.** No part of this manual may be copied, reproduced, translated or reduced to any electronic medium or otherwise without the prior written consent of the copyright holder.

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**EN****Velleman® Service and Quality Warranty**

Since its foundation in 1972, Velleman® acquired extensive experience in the electronics world and currently distributes its products in over 85 countries.

All our products fulfil strict quality requirements and legal stipulations in the EU. In order to ensure the quality, our products regularly go through an extra quality check, both by an internal quality department and by specialized external organisations. If, all precautionary measures notwithstanding, problems should occur, please make appeal to our warranty (see guarantee conditions).

**General Warranty Conditions Concerning Consumer Products (for EU):**

- All consumer products are subject to a 24-month warranty on production flaws and defective material as from the original date of purchase.
- Velleman® can decide to replace an article with an equivalent article, or to refund the retail value totally or partially when the complaint is valid and a free repair or replacement of the article is impossible, or if the expenses are out of proportion. You will be delivered a replacing article or a refund at the value of 100% of the purchase price in case of a flaw occurred in the first year after the date of purchase and delivery, or a replacing article at 50% of the purchase price or a refund at the value of 50% of the retail value in case of a flaw occurred in the second year after the date of purchase and delivery.

**• Not covered by warranty:**

- all direct or indirect damage caused after delivery to the article (e.g. by oxidation, shocks, falls, dust, dirt, humidity...), and by the article, as well as its contents (e.g. data loss), compensation for loss of profits;
- consumable goods, parts or accessories that are subject to an aging process during normal use, such as batteries (rechargeable, non-rechargeable, built-in or replaceable), lamps, rubber parts, drive belts... (unlimited list);
- flaws resulting from fire, water damage, lightning, accident, natural disaster, etc....;
- flaws caused deliberately, negligently or resulting from improper handling, negligent maintenance, abusive use or use contrary to the manufacturer's instructions;
- damage caused by a commercial, professional or collective use of the article (the warranty validity will be reduced to six (6) months when the article is used professionally);
- damage resulting from an inappropriate packing and shipping of the article;
- all damage caused by modification, repair or alteration performed by a third party without written permission by Velleman®.
- Articles to be repaired must be delivered to your Velleman® dealer, solidly packed (preferably in the original packaging), and be completed with the original receipt of purchase and a clear flaw description.
- Hint: In order to save on cost and time, please reread the manual and check if the flaw is caused by obvious causes prior to presenting the article for repair. Note that returning a non-defective article can also involve handling costs.
- Repairs occurring after warranty expiration are subject to shipping costs.
- The above conditions are without prejudice to all commercial warranties.

**The above enumeration is subject to modification according to the article (see article's manual).**

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